

Product description

The NCI/ADR-RES cell line is a multidrug-resistant human ovarian cancer cell line derived from the OVCAR-8 parental line. It is widely used in cancer research to investigate mechanisms of chemoresistance, particularly those involving the overexpression of the MDR1 gene, which encodes the drug efflux transporter P-glycoprotein. This cell line is part of the NCI-60 panel developed by the U.S. National Cancer Institute and serves as a valuable model for evaluating anticancer agents and studying strategies to overcome multidrug resistance. Notably, NCI/ADR-RES was previously misidentified as a derivative of the MCF-7 breast cancer cell line and referred to as MCF-7/ADR, a misclassification that was later corrected through molecular and genomic profiling.

Name: NCI-ADR/RES cell line

Organism: Human

Disease: Cancer

Cancer detailed: High grade ovarian serous adenocarcinoma

Production detail: The MCF-7/ADR cell line was first described in 1986 by Batist et al., who developed it by exposing the MCF-7 breast cancer cell line (mis-identified) to increasing concentrations of the chemotherapeutic drug adriamycin (doxorubicin). This process led to the selection of a stable, multidrug-resistant subline that overexpressed P-glycoprotein (MDR1/ABCB1), making it a widely used model for studying drug resistance in cancer.

Tissue: Ovary

Donor: Female, 64 Years

Parent cell line: OVCAR-8

Growth properties: Adherent

Cellosaurus ID: (CVCL_1452)

Biosafety level: 1

Contributor(s)

Inventor: Kenneth H. Cowan

Institute: National Cancer Institute

Properties

Product format: Frozen

Unpacking and storage:

1. Check all containers for leakage or breakage.
2. Remove the frozen cells from the dry ice packaging and immediately place the cells at a temperature below -130°C, preferably in liquid nitrogen vapor, until ready for use.

Recommended medium: RPMI 1640 + 10% FBS + 50 U/mL Penicillin + 50 µg/mL Streptomycin + 1.84 µM Doxorubicin (to maintain resistance phenotype).

Culture conditions: 37.0°C ± 1.0°C humidified incubator with 5.0% CO₂

Cryopreservation medium: 10% DMSO in FBS

Handling instructions

1. Please ensure that vials are frozen when received, and store at **<-130 °C long term**. When removing frozen cells from storage, it is important to minimize exposure to room temperature (15 - 25°C). If not proceeding directly to thawing, place the cells on dry ice or in a liquid nitrogen container.
2. **Do not thaw at room temperature.** To thaw, swirl the vial quickly in a 37 °C water bath with O-ring and cap above the water to avoid contamination. Remove from the water bath with a small ice pellet remaining (this should not take more than 2 minutes) and wipe the exterior with 70% ethanol or isopropanol before transferring to a biosafety cabinet. Further steps should be conducted under aseptic conditions.
3. We strongly recommend that the volume of cell suspension is measured, and a 20 µL aliquot be set aside at this point for a viable cell count using trypan blue or similar dye.
9. Transfer contents to a 15 mL conical tube containing 9 mL of warm complete medium.
10. Centrifuge at 1000 rpm for 5 minutes to remove DMSO.
11. Discard supernatant and gently resuspend the pellet in fresh complete medium.
12. Seed cells into a T25 or T75 flask with 5–10 mL of medium. Place in 37°C, 5% CO₂ incubator. Change medium after 24 hours to remove residual DMSO and dead cells.
13. Subculture routine: Split 1:3 to 1:6 at ~80% confluency (every 2-3 days) using 0.25% Trypsin-EDTA for detachment at 37 °C for 5 minutes.

References

- Kunkel, et al. Cancer Res. 2024 Aug 1;84(15):2403-2416. PMID: 38861359
- Gholami, et al. Cell Rep. 2013 Aug 15;4(3):609-20. PMID: 23933261
- Ke, et al. Med Oncol. 2011 Dec;28 Suppl 1: S135-41. PMID: 21116879
- Liscovitch and Ravid. Cancer Lett. 2007 Jan 8;245(1-2):350-2. PMID: 16504380
- Adams, et al. J Transl Med. 2005 Mar 4;3(1):11. PMID: 15748285
- D A Scudiero, et al. J Natl Cancer Inst. 1998 Jun 3;90(11):862. PMID: 9625176
- O'Connor PM, et al. Cancer Res. 1997 Oct 1;57(19):4285-90. PMID: 9331090
- Batist G, et al. J Biol Chem. 1986 Nov 25;261(33):15544-9. PMID: 3782078

Material Citation

If use of this material results in a scientific publication, please cite the material in the following manner:
NCI/ADR-RES cell line, was invented by Kenneth H. Cowan (CancerTools.org #161022).

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

While CancerTools.org has made all reasonable efforts to ensure that the information provided by CancerTools.org and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.